

Safety Considerations

2-925-01

Instruction manual

Mini Test Tube Mixer

MVM -10



Thank you for your purchase.

Thank you for your selecting Mini Test Tube Mixer MVM-10.





For your safety & life span of the product, please read this "instruction manual" fully before using and keep these instructions.

After reading this instruction, keep it in a place where you can see it at any time. Make sure that this instruction with the product should be given to the following users for safe use.

AS ONE Co., Ltd.

Please read "Safety considerations" before using and use the product correctly. The suggestions below have possible to connect the serious result.
Please keep these instructions for safety.

Warning mark & meaning

 Danger	Danger of death or serious injury is near at hand with wronghandling.
 Warning	Riskiness of death or serious injury may occur with wronghandling.
 Caution	Injury and physical loss only may occur with wrong handling.
 Suggestion	Caution is required for safety.

Danger

- Do not disassemble or modify the equipment. It may cause electric shock or accident.
- Since this product operates at high speed, be careful not to let your hair or other objects get caught inside. Also, please do not touch the pad directly as it is dangerous.


Caution

- Be sure to hold the plug firmly when pulling out the power cord. Also when inserting the plug, make sure that it is plugged in.
 - This product is designed for indoor use. Do not use outdoors or in areas exposed to water.
 - Keep the ambient temperature within the range of 5 ~ 35°C
- Also, to avoid malfunction, do not use this product in direct sunlight or near heating appliances .
- please be fully careful of power supply voltage.

•Specifications

Model	MVM -10
Rotation speed	800 ~ 2500 rpm
Operating mode	Continuous Operation Type/Touch Start Type
Size/Weight	72 x 88 x 82 mm/Approximately 0.3 kg
Power supply	AC100V to 240V 50/60 Hz (dedicated AC adapter) Or AA size battery x 4 pieces (with battery case sold separately)
Accessories	Dedicated AC adapter

•Name of parts

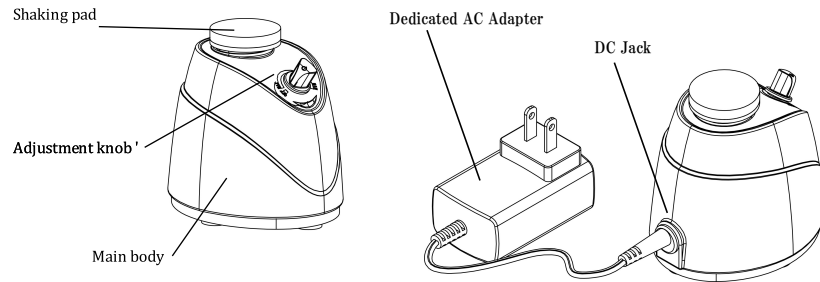

•How to use

- ① Place the product on an empty horizontal position within 10cm of the circumference.
- ② Set the adjustment knob to "OFF", and connect the AC adapter to the power supply.

•Continuous Operation mode

When the adjustment knob is rotated in clockwise, the shaking pad will shake. Please adjust the rotation speed according to the MAX-MIN display, and stir the sample.

•Touch Start Mode

If the shaking pad is pressed from the top when the adjustment knob is set to [OFF], operation starts in "touch start mode".

In this mode, the rotation speed will be fixed (Approximately 2500 rpm).

•The machine can also be operated in battery using an optional battery case (2-925-11)
Also, please replace the battery when the rotation speed is low during use.

•Product Maintenance

- Do not use this product in the presence of corrosive gas or liquid. Do not use organic solvents for maintenance, and use a soft cloth when you need to wipe off dirt.
- If there was an abnormal noise, odor, or other abnormalities occur during use, discontinue using this product immediately and contact your dealer.
- Store this product in a well-ventilated area.
Storing in a humid environment can cause damage.
- When the shaking pad is worn, it can be replaced. Purchase an optional replacement attachment hammer (Ason Code: 2-925-12).
(This pad can be adhered with the double-sided tape inside.)

For further information, please contact:

Department in charge : AS ONE Corporation Overseas Sales Department

Address : 2-1-27,Edobori,Nishi-ku Osaka550-8527 Japan

TEL:81-6-6447-8905

FAX:81-6-6447-8681

E-mail: asoneint@so.as-1.co.jp

Made in China
March 2021 Published edition
of J2_ (1)